

# Assembly Instructions: Fiberglass Back Mount Lab Frame Kit, 8 Foot

## Lattice Frame:

Use the closed connectors and lattice rods to assemble the frame. Use one open connector at each rod intersection. The kit assumes a lattice spacing of 1-foot; however, this may vary depending on your application. Rods may be cut for fit as necessary.

## Mounting:

After assembling the lattice frame, there should be an equal number of remaining 4-inch rods, baseplates, and open connectors. Connect an open connector to one end of the 4-inch rod and a base plate to the other end. Attach the open connector to the lattice frame 90 degrees in a square or rectangular pattern (the base plate pointing to the back mounting surface). Base plates may be secured with #10 screws. Mounting Screws Not Included.

*NOTE: Follow your fume hood manufacturers' requirements for attachment.*

*\*Images may be different than actual product*



36" Rod X 7

86" Rod X 2

4" Rod X 8



X 22



X 8

