

# Assembly Instructions: Aluminum Table Top Lab Frame Kit, 8 Foot

## Lattice Frame:

Use the connectors and lattice rods to assemble the frame. The kit assumes a lattice spacing of 1-foot and a border on the sides and top. Spacing may vary depending on your application. Rods may be cut to fit as necessary.

## Mounting:

After assembling the lattice frame attach vertical base plates to the bottom of the vertical rods. There should be an equal number of remaining 24" rods, swivel baseplates and connectors. Connect a connector to one end of the 24" rod and a swivel base plate to the other end. Attach the connector to the lattice frame 30-45 degrees. Base plates may be secured with #10 screws. **Mounting Screws Not Included.**

*NOTE: Follow your fume hood manufacturers' requirements for attachment.*

*\*Images may be different than actual product*




36" Rod X 8

86" Rod X 3

24" Rod X 4



X 28

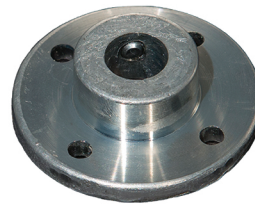

X 8



X 4

